



## Atomx Firmware - 102.1298.atomxmain.S19

---

This latest firmware upgrade includes enhancements to increase overall functionality of the unit: (see readme file included in 102.1298.atomxmain.S19.zip below for more details)

### Installation Instructions (firmware) ▼

---

- 1.) Right click the Download Now (Firmware) link.
- 2.) Select "Save Target As..."
- 3.) In the "Save As" prompt, you should see the name of the file (102.1298.atomxmain.s19) in the "File Name" text box. In the dropdown menu of "Save as type," select "All files."
- 4.) Save the file (102.1298.atomxmain.s19) in the default directory (C:\Program Files\Teledyne Tekmar\Atomx TekLink\Firmware Updates).
- 5.) Open VOC TekLink and flash the unit while running TekLink.
- 6.) For more information on "flashing", go to TekLink Help, type in "Flash" and follow the links for flashing.